

INSTRUCTIONS FOR REPLACING

ViewPoint™

By MISCO®

BRINGING REFRACTOMETERS
OUT OF THE DARK AGES...

The ViewPoint Illuminator provides light for the refractometer, so the scale can be read in dim light or total darkness. With the ViewPoint, you will not need to point the refractometer at another light source.

1. Remove Hinge Pin

Remove the hinge pin by carefully pushing it part way through the cover plate with a small nail or pinch. When enough of the pin is exposed on one side of the cover plate, hold the cover plate to the body of the refractometer with one hand, while gripping the exposed pin using pliers in your other hand. Pull the pin the rest of the way out. Be careful not to bend the pin. Also, be careful not to lose the small ball bearing and spring located under the left leg of the ViewPoint.



2. Replace Cover Plate

With the ball bearing/spring assembly in place, reinsert the pin through the body and ViewPoint by reversing the steps in point 1 above.

For Technical Assistance, Please Call:
800-358-1100